OmniCure®

X-Cite®

Safety Precautions

035-00636R Rev 0

[ES] Medidas de seguridad.....	1
[CS] Bezpečnostní opatření.....	8
[DA] Sikkerhedsforanstaltninger.....	15
[DE] Sicherheitsvorkehrungen.....	22
[EN] Safety Precautions	29
[FR] Consignes de sécurité	35
[IT] Precauzioni di sicurezza.....	42
[NL] Veiligheidsmaatregelen.....	49
[NO] Sikkerhetsforanstaltninger.....	56
[PL] Środki bezpieczeństwa.....	62
[PT] Precauções de Segurança.....	69
[FI] Turvaohjeet.....	76
[SV] Säkerhetsföreskrifter.....	82

[ES] Medidas de seguridad

Glosario de símbolos

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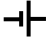
[EN] Safety Precautions

Glossary of Symbols

Symbol	Meaning	Symbol	Meaning
	CAUTION - if the precaution is not taken, it may cause minor or moderate injury.		BATTERY
	WARNING – if the warning is not heeded, it can cause death or serious injury.		
	DANGER – if the danger is not avoided, it will cause death or serious injury.		

General Safety Precautions

CAUTION

- (*X-Cite only*) Use of the X-Cite system is intended only for microscopes using Excelitas Technologies collimating adaptors. Any attempt to use this system with a microscope not approved by Excelitas Technologies is not recommended and may impair system safety.
- To prevent damage/degradation of the light guide, allow adequate clearance at the front/rear of the unit to prevent kinking or excessive bending.
- If this unit is used in a manner not specified by Excelitas Technologies, the protection provided by the equipment may be impaired.
- Any electronic equipment connected to this product must comply with the requirements of EN/IEC 60950.

WARNING

- UV emitted from this product. Avoid eye and skin exposure to unshielded product. Do not look at operating lamp/LED. Eye injury may result. Always use protective eyewear and turn the unit off before removing any cover.
- Never look into the light emitting end of the light guide/LED head. The light could severely damage the cornea and retina of the eye if the light is observed directly. Eye shielding must be used at all times as well as clothing to protect exposed skin.

- Make sure the light guide/LED head is securely attached to the proper mounting fixture, as described in the User Guide, prior to turning on power to the unit. This will minimize the risk of exposure to the UV light.
- Only use this product with the external power supply (if applicable) and power cord set provided. Ensure the external power supply cord is connected to a grounded 3-pin outlet only. Any substitution of these components will invalidate regulatory certification of this product and may impair operating safety.
- To reduce the risk of fire or shock, always replace the fuses with the same type and rating. Always detach the power supply cord prior to attempting to replace fuses.
- To disconnect the main supply source, put the on/off button in the off position and unplug the power cord.
- This unit is designed for bench top use only. Ensure that the unit is placed on a hard, stable surface, ensuring that ventilation openings are not obstructed. Any obstruction of these openings could result in a possible over-heating condition.
- Manual operation:
The level of UV and visible energy supplied by this product is sufficient to ignite flammable substances. During manual operation, the unit must be attended at all times by a qualified operator. The unit must not be left unattended while turned on. If an operator leaves the work area of the unit, the power switch must be turned off.
- Automated operation:
The level of UV and visible energy supplied by this product is sufficient to ignite flammable substances. When the unit is operated unattended in an automated environment, an alarm function must be provided by the user to indicate a malfunction in the associated equipment used.

OmniCure and X-Cite Lamp-Based Systems

CAUTION

- The lamp module's operational life can be significantly shortened if it is handled incorrectly. Do not touch the bulb's glass envelope or the inside surface of the reflector. Skin oils can cause the lamp module to fail prematurely.
- Allow the lamp module to cool down completely before opening the unit and handling the lamp module.

 **WARNING**

- IR emitted from this product. Do not look at operating lamp.
- Hg – LAMP CONTAINS MERCURY, manage in accordance with disposal laws, see: www.lamprecycle.org or 1-800-668-8752.

DANGER: Exposure to Mercury represents a health hazard to humans.

When unpacking or installing the lamp, always wear protective clothing and a face mask. Operate the lamp inside the unit's lamp housing only. This prevents direct viewing of the arc and, in the case of lamp bursting, contains the lamp particles. In the rare instance in which a lamp bursts and the mercury content is released, the following safety precautions are recommended: all personnel should be immediately evacuated from the area to prevent inhalation of the mercury vapor. The area should be well ventilated for a minimum of 30 minutes. Prior to clean up ensure an "approved-for-mercury" respirator mask and non-porous gloves such as latex or rubber are used. After the lamp housing elements have cooled, the mercury residue should be collected with the use of a special absorbing agent available from laboratory equipment suppliers. Listed below are examples of web sites for obtaining Mercury Spill Kits:

<http://rosshc.org/mercury-safety/>

http://www.coleparmer.ca/Category/MERCSORB_Mercury_Spill_Kit/43577

- The method in which lamps are disposed of must comply with local rules & regulations for disposal of hazardous materials. Lamps may be returned to Excelitas Technologies providing they are returned in their original packaging. Excelitas Technologies will dispose of them in the appropriate manner.

 **DANGER**

- This unit contains HIGH VOLTAGE components. It is recommended that ONLY QUALIFIED TECHNICAL PERSONNEL perform any testing or repairs of this unit. If servicing the unit, disconnect the external power supply before opening the cover of this unit. All cover screws must be replaced prior to applying power to the unit or safety of the unit will be impaired. Note that opening the cover and/or altering the unit in any way may void the warranty.

OmniCure LED-Based Systems

CAUTION

- Do not attempt to connect any other type of equipment to the connectors labelled “CH1-2” and “CH1-4” located on the rear panel of the OmniCure LX180 and LX400+, respectively, or any other multi-head unit. These connectors are provided only for connection of the various LED heads available from Excelitas Technologies.
- Due to elevated operating temperatures, avoid contact of the LED head(s) when energized (front panel “LED ON” emission indicator is illuminated red). The LED heads are designed to be mounted in a suitable fixture prior to use. User clamp type heat sink assemblies are provided with each LED head to provide user safety and optimum thermal management. Prior to handling of the UV LED head(s), allow a cool down for a period of approximately five minutes after system power has been removed.
- The UV LED head’s operation can be affected if handled incorrectly. Never touch the protective lens cover assembly of the UV LED head. The presence of skin oils may result in a decrease in system performance.
- To protect the system from electrostatic-related damage, which can disable the system or cause permanent damage or require repair by service technicians, place all modules on an antistatic surface, such as an antistatic discharge mat, an antistatic bag, a disposable antistatic mat or a grounded metal surface. If possible, wear an antistatic grounding strap connected to a metal surface on the chassis when working on the system.

WARNING

- The protective eyewear provided is designed only for viewing of indirect UV light on an intermittent basis. Consistent with this intended use, this eyewear meets the exposure limits as defined by HSM-73-11009 from the U.S. Department of Health, National Institute for Occupational Safety and Health. Additional protective UV shielding is recommended for any process that requires direct viewing or visual monitoring of UV light for extended periods of time. It is the responsibility of the user to ensure that adequate safeguards are implemented to prevent any potential hazardous UV exposure.
- The LEDs available for this system provide optical output power classified as RISK GROUP 2 per IEC/ EN 62471-2. The user-related risk of this system depends on the final installation and use of this product as detailed in the User Guide.

- Handheld use of the LED heads is not recommended and may expose the user to dangerous optical radiation. Additionally, to prevent unintentional exposure, use of the on/off key switch is recommended.
- Ensure the system controller power is turned off before disconnecting or re-connecting any of the UV LED heads.
- Wear/use proper UV shielding/protection within 46cm of the UV LED heads (if applicable).

OmniCure R2000 Radiometer

X-Cite Optical Power Measurement System (OPMS: XR2100 & XP750)

CAUTION

- In instances where high power light sources are measured for extended periods of time, the light guide adaptors supplied with the R2000/XR2100 may become hot. Use caution when handling these adaptors.
- Allow adequate clearance to prevent kinking or excessive bending to prevent damage/degradation of the attached sensor cable.
- The detection window operational life can be significantly shortened if handled incorrectly. Do not touch the window surface. Skin oils and scratches can cause the window to fail prematurely.
- Do not discard used batteries. Return to the nearest authorized Excelitas Technologies service center for disposal/recycling.

WARNING

- Make sure the light guide is properly inserted into the R2000/XR2100 and/or the microscope (if applicable) before working with the system. This will minimize the risk of exposure to the UV light.
- The R2000/XR2100 is supplied with two lithium batteries. Lithium batteries present a potential fire, explosion or severe burn hazard. DO NOT attempt to re-charge, disassemble, incinerate, short circuit or expose battery to temperatures above 100°C or expose contents to water.

BATTERY

- Lithium batteries must have terminals taped with non-conductive material prior to returning for disposal/recycling to prevent short-circuiting. External packaging material must provide adequate protection to contents.

- The lithium batteries supplied in R2000/XR2100 DO NOT contain: mercury, lead, manganese or cadmium. Substitution of any other type of battery is not recommended and may void warranty.

X-Cite XLED1

CAUTION

- Do not attempt to connect the Touch Screen Controller to any type of electrical equipment other than the XLED1 main system.
- To protect the LED module, driver and system from electrostatic-related damage, which can disable the system, permanently damage the LED module and driver, or require repair by service technicians, place all modules on an antistatic surface, such as an antistatic discharge mat, an antistatic bag, a disposable antistatic mat or a grounded metal surface. If possible, wear an antistatic grounding strap connected to a metal surface on the chassis when you replace any LED module or LED driver.

WARNING

- The service cover should be removed by qualified technical personnel only. No user serviceable parts are located within the XLED1 chassis or the Touch Screen Controller. This system contains hazardous voltages—do not attempt to remove access covers while equipment is energized.
- Do not lift the XLED1 from the rear of the removable housing when system is either on or off.

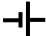
X-Cite TURBO

WARNING

- The X-Cite TURBO is classified as Class 1 laser product. IEC 60825-1: Safety of laser products – Part 1: Equipment classification and requirements: 2014-05. This class is eye-safe under all operating conditions. This class includes all lasers or laser systems which cannot emit levels of optical radiation above the exposure limits for the eye under any exposure conditions inherent in the design of the laser product. There may be a more hazardous laser embedded in the enclosure of a Class 1 product, but no harmful radiation can escape the enclosure.

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Excelitas Canada Inc.